



SAFETY ENCLOSURES, CLEAN BENCHES &
NON-VENTILATED ENCLOSURES






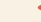


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
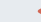
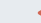
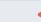






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
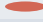

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


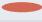
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


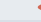
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
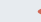



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
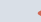
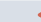
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



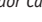
















Protecting your
laboratory environment

LABCONCO

SAFETY ENCLOSURES & CLEAN BENCHES

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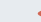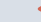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SELECTION GUIDE

Description Product	Class I Purifier Class I, Safety Enclosure See pages 2-7	 Purifier HEPA Filtered Enclosure See pages 2-7	Class I with Purifier Class I, Safety Enclosure See pages 2-7	 Canopy Duct Purifier HEPA Filtered Enclosure See pages 2-7
Protection Provided Product Only Personnel and Environment				
Suitable For Use With: Biosafety Level 1, 2 and 3 Agents Objectionable Odors Toxic Particulates Volatile Toxic Chemicals Tracer Quantities of Radionuclides	  	  	   	   
Airflows & Exhaust System Exhausts to: Exhaust Duct Connection Ganged Exhaust System Possible % Recirculation In Cabinet	ROOM NONE 	ROOM NONE 	OUTSIDE HARD DUCT  	OUTSIDE HARD DUCT  
Options/Accessories Nominal Width Service Fixtures UV Light IV Bar Work Surface Telescoping or Hydraulic Lift Base Stand Vibration Isolation Table Ductwork/Remote Blower Trace Odor Carbon Filters	2', 3', 4'     	2', 3', 4'    	2', 3', 4'     	2', 3', 4'    
APPLICATIONS				
Biotechnology Sterile Media Preparation Non-biohazardous Culture Maintenance Non-biohazardous Tissue Culture Maintenance Plant Tissue Culture Maintenance Blood Component Research Human Tissue Research PCR				
Microbiology Sterile Media Preparation Non-biohazardous Culture Maintenance Culture Maintenance Work with Malodorous Cultures Non-biohazardous Tissue Culture Maintenance Non-volatile, Non-toxic Staining Staining with Volatile, Toxic Material Non-volatile Radiolabeling/Tagging Volatile Radiolabeling/Tagging	  	  	   	   
Pharmacy Non-toxic IV Solution Preparation				
General Tissue Fixation/Staining Preparation Weighing Toxic Powders Asbestos Handling Suspicious Mail Handling Electronics Inspection/Repair Soldering	    	    	   	   

*When Trace Odor Carbon Filter is used

SELECTION GUIDE

[illegible]

PURIFIER® CLASS I & HEPA FILTERED SAFETY ENCLOSURES

I N T R O D U C T I O N

Purifier Class I Safety Enclosures and Purifier HEPA Filtered Enclosures protect you and your laboratory environment. They offer an economical alternative to Class II laminar flow cabinets when your applications do not require product protection.

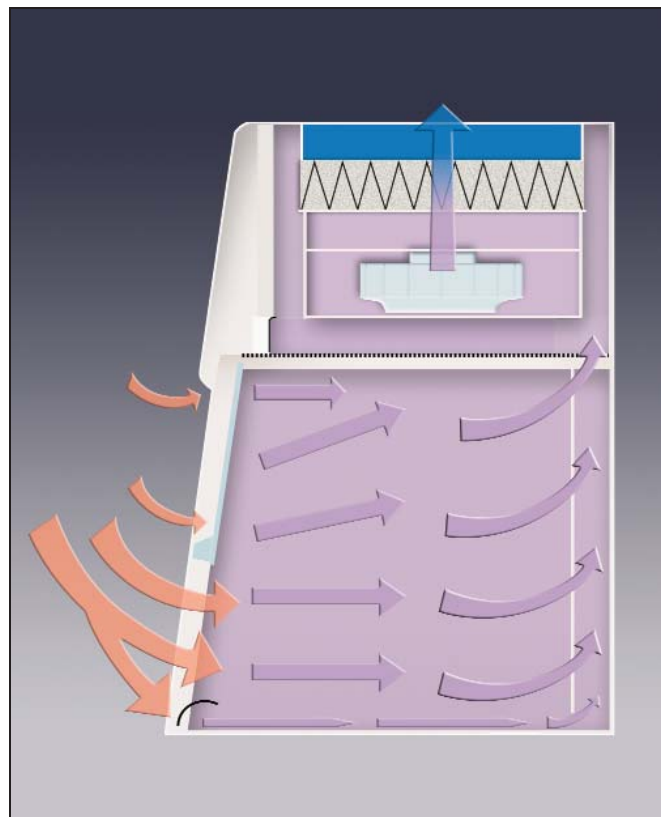
Purifier Class I Enclosures protect the user from agents that require Biosafety Level 1, 2 or 3 containment but no product protection. These enclosures include an ultraviolet light that may be used in conjunction with surface disinfection to ensure thorough decontamination. The UV light is turned on only while the enclosure is unattended, to avoid personnel exposure to UV rays.

Purifier HEPA Filtered Enclosures, which do not include a UV light, protect the user from hazardous chemical powders, dust and allergens in diverse applications ranging from screening suspicious mail to blending dry chemicals.


For applications involving odor-generating materials, Purifier Class I and HEPA Filtered Enclosures may be outfitted with an accessory Trace Odor Carbon Filter to trap nuisance organic, formaldehyde or ammonia vapors. When connected to the accessory Canopy Connection and ducted outside to a remote blower or house exhaust, Purifier Class I and HEPA Filtered Enclosures may be used with fume-generating materials or agents that have been treated with toxic chemicals or radionuclides as an adjunct to research. These canopy-ducted enclosures have the added benefit of eliminating nuisance odors. Use of acids, which would corrode the internal blower within these enclosures, is prohibited. Contact Labconco for information on enclosures suitable for use with acids.



Purifier HEPA Filtered Enclosure 3980302 on Solid Epoxy Dished Work Surface 3909901 and Telescoping Base Stand 3746701 with Ergonomic Chair 3744000.



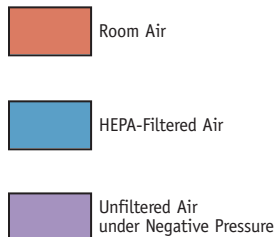During operation, room air is drawn into the front of the enclosure, preventing aerosols from escaping into the room. Before leaving the enclosure and returning to the environment, 100% of the air passes through an exhaust HEPA filter.




PURIFIER® CLASS I & HEPA FILTERED SAFETY ENCLOSURES


FEATURES & BENEFITS


Sturdy epoxy-coated aluminum frame and steel rear plenum and baffle dissipate static charge and provide durability and corrosion resistance.

Front-mounted Minihelic pressure gauge monitors pressure differential across the HEPA filter to help determine filter loading.

Tempered safety glass front sash and sides offer excellent visibility and protection. Glass provides better fire, scratch and corrosion resistance than acrylic and dissipates static charge.

 **Upper containment sash foil** bleeds air into the enclosure to direct contaminants away from the operator's breathing zone.

 **Patented* rear baffle with zones of perforations** promotes horizontal laminar airflow to maximize containment.

 **Side-entry air foils** allow air to sweep across the interior glass surfaces to enhance containment.

Two utility ports with iris openings allow passage of tubing, electrical cords and cables from inside to back exterior for connection to services.

Accessory Telescoping Base Stand. See pages 24-25 for ordering information.

Particulate containment tested. Free validation package containing independent containment testing verification is available.

ASHRAE 110-1995 conformity.

Space-saving design. Impeller and HEPA filter are contained within the enclosure's upper plenum for a smaller overall footprint than a separate enclosure and exhaust can offer.


Built-in exhaust HEPA filter, 99.99% efficient on 0.3 micron particles.

Intrinsically-safe design. The positive pressure impeller is surrounded by a negative pressure plenum so that if a leak should develop, the unfiltered air is captured and refiltered.

Optional built-in Guardian™ Airflow Monitor alerts the user to face velocities outside of preset range, indicating HEPA filter loading or blower malfunction.

Glare-free fluorescent lighting.

Low discharge volume reduces cross currents and provides quiet operation from 49 to 67 decibels depending on face velocity.

 **Upper dilution air supply** introduces room air at the top of the sash to dilute concentrations in the upper chamber and bathes the back of the sash with clean air.

Front-mounted light and blower switches.

Speed control regulates speed of the built-in blower and may be used by a certifier to adjust and validate inflow velocity.

Ergonomic 10° angled sash allows a closer view, reduces glare and provides a more comfortable operating position than vertical sashes. The hinged sash pivots upward and locks to a loading height of 20". Side seals help to keep contaminants contained.

 **Patented* Clean-Sweep™ air foil** allows air to sweep the work surface for maximum containment. Airflow openings pull inflow air from under the air foil so that clean air continually flows over the air foil creating a constant barrier of protection from contaminants.

Accessory Solid Epoxy Dished Work Surface. See page 6.

ETL-listed. Enclosures carry the ETL mark signifying that they are certified to UL 3101-1/61010-1 and CAN/CSA C22.2 No. 1010.1.

CE mark. All 230 volt, 50 Hz models conform to the CE (European Community) requirements for electrical safety and electromagnetic compatibility.

Optional Trace Odor Carbon Filters designed for organic, formaldehyde and ammonia odor applications are easily installed and replaced from the top of the enclosure.



Biological containment verified per NSF Standard Number 49 Personnel Protection Test.

254 nm UV light (Class I models only) for secondary decontamination while the enclosure is not in use.



* U.S. Patent No. 6,461,233

 Exclusive feature

PURIFIER® CLASS I SAFETY ENCLOSURES

ORDERING INFORMATION

Standard Features

- Nominal inflow velocity of 75 to 105 fpm
- Ergonomic air foil with aerodynamic Clean-Sweep* airflow openings
- Upper dilution air supply
- Zone-perforated rear baffle*
- Side-entry air foils
 - Exhaust HEPA filter, 99.99% efficient on 0.3 micron particles
 - Variable speed built-in blower with dynamically-balanced, vibration-isolated motorized impeller and solid state control rated for 5 amps
- Intrinsically-safe, negative pressure design
- Minihelic pressure gauge
- Low discharge volume and quiet operation (*see chart below*).

Nominal Width	Total Exhaust CFM and Noise Pressure					
	105 fpm	db(A)	90 fpm	db(A)	75 fpm	db(A)
2 Feet	155	54-59	130	51-57	110	49-55
3 Feet	230	60-64	200	55-58	165	52-56
4 Feet	310	64-67	265	60-63	220	54-58

- Upper containment sash foil
 - Glacier white and gray epoxy-coated aluminum frame and steel rear plenum and baffle
 - Fluorescent lamp
 - 254 nm UV lamp
 - Blower switch and 3-way light switch: UV, off and fluorescent
 - 1/4" thick tempered safety glass front sash and sides
 - 9.44" high sash opening
 - Ergonomic 10° angled and hinged sash that pivots upward, locks to a loading height of 20" and has side wiping seals to contain contaminants
- Two utility ports with iris openings
- Particulate containment tested
- Biological containment verified per NSF Standard Number 49 Personnel Protection Test
- ASHRAE 110-1995 conformity tested
- ETL listing, UL 3101-1/61010-1 and CAN/CSA C22.2 No. 1010.1 conformity (115 volt models)
- CE Conformity Marking (230 volt models)
- One year warranty on parts and service
- Validation package available



Purifier Class I Safety Enclosure 3980203 on Solid Epoxy Dished Work Surface 3909900 and Telescoping Base Stand 3746700.

Standard Option

- Guardian Airflow Monitor that continuously monitors airflow and has two LED lights to indicate when airflow is within or outside set point range. The green LED glows when airflow is within set point range. The red LED glows and an audible alarm sounds when airflow is outside set point range.

Required Accessories

- Supporting work surface. Place on existing casework or see page 6.
- Supporting base. Place on existing casework or see pages 24-27.

See page 6 for dimensional drawings.

Catalog #	Nominal Width	Exhaust Volume	Electrical Requirements	Airflow Monitor	Power Cord & Plug	Shipping Weight
3980201	2 feet	110 - 155 CFM	115 volts, 60 Hz, 3 amps		115 volts, 15 amps	125 lbs. (57 kg)
3980203	2 feet	110 - 155 CFM	115 volts, 60 Hz, 3 amps	•	115 volts, 15 amps	125 lbs. (57 kg)
3980221†	2 feet	110 - 155 CFM	230 volts, 50 Hz, 2 amps		230 volts, 10 amps, no plug	125 lbs. (57 kg)
3980223†	2 feet	110 - 155 CFM	230 volts, 50 Hz, 2 amps	•	230 volts, 10 amps, no plug	125 lbs. (57 kg)
3980301	3 feet	165 - 230 CFM	115 volts, 60 Hz, 3 amps		115 volts, 15 amps	160 lbs. (73 kg)
3980303	3 feet	165 - 230 CFM	115 volts, 60 Hz, 3 amps	•	115 volts, 15 amps	160 lbs. (73 kg)
3980321†	3 feet	165 - 230 CFM	230 volts, 50 Hz, 2 amps		230 volts, 10 amps, no plug	160 lbs. (73 kg)
3980323†	3 feet	165 - 230 CFM	230 volts, 50 Hz, 2 amps	•	230 volts, 10 amps, no plug	160 lbs. (73 kg)
3980401	4 feet	220 - 310 CFM	115 volts, 60 Hz, 3 amps		115 volts, 15 amps	195 lbs. (88 kg)
3980403	4 feet	220 - 310 CFM	115 volts, 60 Hz, 3 amps	•	115 volts, 15 amps	195 lbs. (88 kg)
3980421†	4 feet	220 - 310 CFM	230 volts, 50 Hz, 2 amps		230 volts, 10 amps, no plug	195 lbs. (88 kg)
3980423†	4 feet	220 - 310 CFM	230 volts, 50 Hz, 2 amps	•	230 volts, 10 amps, no plug	195 lbs. (88 kg)

• Exclusive feature

* U.S. Patent No. 6,461,233

† International electrical configuration

PURIFIER® HEPA FILTERED ENCLOSURES

ORDERING INFORMATION

Standard Features

- Nominal inflow velocity of 75 to 105 fpm
- Ergonomic air foil with aerodynamic Clean-Sweep* airflow openings
- Upper dilution air supply
- Zone-perforated rear baffle*
- Side-entry air foils
- Exhaust HEPA filter, 99.99% efficient on 0.3 micron particles
- Variable speed built-in blower with dynamically-balanced, vibration-isolated motorized impeller and solid state control rated for 5 amps
- Intrinsically-safe, negative pressure design
- Minihelic pressure gauge
- Low discharge volume and quiet operation (*see chart below*).

Nominal Width	Total Exhaust CFM and Noise Pressure					
	105 fpm	db(A)	90 fpm	db(A)	75 fpm	db(A)
2 Feet	155	54-59	130	51-57	110	49-55
3 Feet	230	60-64	200	55-58	165	52-56
4 Feet	310	64-67	265	60-63	220	54-58

- Upper containment sash foil
- Glacier white and gray epoxy-coated aluminum frame and steel rear plenum and baffle
- Fluorescent lamp
- Switches for blower and light
- 1/4" thick tempered safety glass front sash and sides
- 9.44" high sash opening
- Ergonomic 10° angled and hinged sash that pivots upward, locks to a loading height of 20" and has side wiping seals to contain contaminants
- Two utility ports with iris openings
- Particulate containment tested
- Biological containment verified per NSF Standard Number 49 Personnel Protection Test
- ASHRAE 110-1995 conformity tested
- ETL listing, UL 3101-1/61010-1 and CAN/CSA C22.2 No. 1010.1 conformity (115 volt models)
- CE Conformity Marking (230 volt models)
- One year warranty on parts and service
- Validation package available



Purifier HEPA Filtered Safety Enclosure 3980402 on Solid Epoxy Dished Work Surface 3909902 and Telescoping Base Stand 3746702.

Standard Option

- Guardian Airflow Monitor that continuously monitors airflow and has two LED lights to indicate when airflow is within or outside set point range. The green LED glows when airflow is within set point range. The red LED glows and an audible alarm sounds when airflow is outside set point range.

Required Accessories

- Supporting work surface. Place on existing casework or see page 6.
- Supporting base. Place on existing casework or see pages 24-27.

See page 6 for dimensional drawings.

Catalog #	Nominal Width	Exhaust Volume	Electrical Requirements	Airflow Monitor	Power Cord & Plug	Shipping Weight
3980200	2 feet	110 - 155 CFM	115 volts, 60 Hz, 3 amps		115 volts, 15 amps	125 lbs. (57 kg)
3980202	2 feet	110 - 155 CFM	115 volts, 60 Hz, 3 amps	•	115 volts, 15 amps	125 lbs. (57 kg)
3980220†	2 feet	110 - 155 CFM	230 volts, 50 Hz, 2 amps		230 volts, 10 amps, no plug	125 lbs. (57 kg)
3980222†	2 feet	110 - 155 CFM	230 volts, 50 Hz, 2 amps	•	230 volts, 10 amps, no plug	125 lbs. (57 kg)
3980300	3 feet	165 - 230 CFM	115 volts, 60 Hz, 3 amps		115 volts, 15 amps	160 lbs. (73 kg)
3980302	3 feet	165 - 230 CFM	115 volts, 60 Hz, 3 amps	•	115 volts, 15 amps	160 lbs. (73 kg)
3980320†	3 feet	165 - 230 CFM	230 volts, 50 Hz, 2 amps		230 volts, 10 amps, no plug	160 lbs. (73 kg)
3980322†	3 feet	165 - 230 CFM	230 volts, 50 Hz, 2 amps	•	230 volts, 10 amps, no plug	160 lbs. (73 kg)
3980400	4 feet	220 - 310 CFM	115 volts, 60 Hz, 3 amps		115 volts, 15 amps	195 lbs. (88 kg)
3980402	4 feet	220 - 310 CFM	115 volts, 60 Hz, 3 amps	•	115 volts, 15 amps	195 lbs. (88 kg)
3980420†	4 feet	220 - 310 CFM	230 volts, 50 Hz, 2 amps		230 volts, 10 amps, no plug	195 lbs. (88 kg)
3980422†	4 feet	220 - 310 CFM	230 volts, 50 Hz, 2 amps	•	230 volts, 10 amps, no plug	195 lbs. (88 kg)

† Exclusive feature



* U.S. Patent No. 6,461,233

† International electrical configuration

PURIFIER® CLASS I & HEPA FILTERED SAFETY ENCLOSURES

D I M E N S I O N S & A C C E S S O R I E S

	2-Foot	3-Foot	4-Foot
A	24.0"	36.0"	48.0"
B	21.6"	33.6"	45.6"

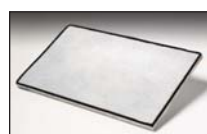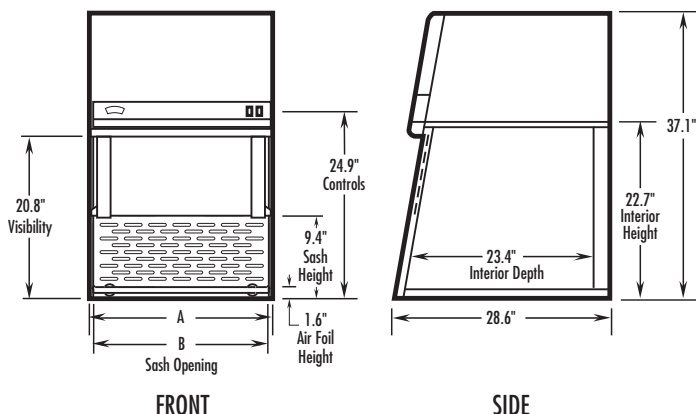

Trace Odor Carbon Filters

For easy installation atop the Purifier Class I or HEPA Filtered Enclosure. Filters trap non-volatile nuisance odors and vapors from samples or minute quantities of chemicals.

Whether or not a Trace Odor Carbon Filter is appropriate for your application requires consideration of factors such as the time-weighted exposure limits of organic solvents, formaldehyde, ammonia or amines. Using Labconco's exclusive Chemical Assessment Program, our technical specialists can help you determine if the Purifier Enclosure with Trace Odor Carbon Filter is right for your applications and, if so, estimate filter life. To begin the process, please complete the Chemical Usage Assessment Form available on our website at: www.labconco.com/chemical_usage.html or call us at 800-821-5525 or 816-333-8811.

Catalog #	Description	For use with:	Shipping Weight
3937200	Organic, 5.5 lbs. activated carbon	2-Ft Purifier Enclosure	6.0 lbs. (2.7 kg)
3937300	Organic, 9.2 lbs. activated carbon	3-Ft Purifier Enclosure	10.0 lbs. (4.5 kg)
3937400	Organic, 12.9 lbs. activated carbon	4-Ft Purifier Enclosure	13.0 lbs. (5.9 kg)
3937201	Formaldehyde, 7 lbs. impregnated carbon	2-Ft Purifier Enclosure	7.5 lbs. (3.4 kg)
3937301	Formaldehyde, 12 lbs. impregnated carbon	3-Ft Purifier Enclosure	12.5 lbs. (5.7 kg)
3937401	Formaldehyde, 16.8 lbs. impregnated carbon	4-Ft Purifier Enclosure	17.0 lbs. (7.7 kg)
3937202	Ammonia & Amines, 8 lbs. impregnated carbon	2-Ft Purifier Enclosure	8.5 lbs. (3.9 kg)
3937302	Ammonia & Amines, 13.5 lbs. impregnated carbon	3-Ft Purifier Enclosure	14.0 lbs. (6.4 kg)
3937402	Ammonia & Amines, 18.9 lbs. impregnated carbon	4-Ft Purifier Enclosure	19.0 lbs. (8.6 kg)



Solid Epoxy Dished Work Surfaces

Black chemical-resistant work surface is contoured to fit the dimensions of the Purifier Class I and HEPA Filtered Enclosure and to contain spills.


Catalog #	For use with	Dimensions	Shipping Weight
3909900	2-Ft Purifier Enclosure	24.0" w x 29.0" d x 1.0" h (61 cm x 73.7 cm x 2.5 cm)	50 lbs. (23 kg)
3909901	3-Ft Purifier Enclosure	36.0" w x 29.0" d x 1.0" h (91.4 cm x 73.7 cm x 2.5 cm)	75 lbs. (34 kg)
3909902	4-Ft Purifier Enclosure	48.0" w x 29.0" d x 1.0" h (122 cm x 73.7 cm x 2.5 cm)	100 lbs. (45 kg)



Base Stands

Epoxy-coated steel. Telescoping Base Stands adjust in 1" increments to provide a working height from 30" to 36" and are available with fixed feet or casters. Electric and Manual Hydraulic Lift Base Stands have infinite

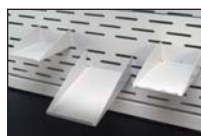height adjustment using a push button or hand crank. See pages 24-27 for more information.



3925000 Utility Shelves

Three epoxy-coated steel shelves attach to the rear baffle perforations to hold a printer, utensils and other miscellaneous supplies. The utensil shelf has a scalloped front edge to support spatulas, brushes and other items. Shipping weight 5.0 lbs. (2.2 kg)



3927700 Bottle Holder

Epoxy-coated steel. Attaches to the rear baffle perforations to hold a squirt bottle or other round container. 3" diameter opening. Shipping weight 1.0 lb. (0.5 kg)



3927800 Tissue Holder

Epoxy-coated steel. Attaches to the rear baffle perforations to hold a 4.63" x 4.75" box of tissue wipers in the upside down position. An opening in the bottom of the holder permits tissues to be dispensed from the bottom. Shipping weight 1 lb. (0.5 kg)



3744000 Ergonomic Chair with Armrests

Chair has 6-way articulating seat and back control for personalized adjustment. Pneumatic mechanism adjusts seat height from 18.25" to 25.75". Shipping weight 35.0 lbs. (15.9 kg). See page 28 for more information.



3746000 Adjustable Footrest

Elevates feet and permits angle repositioning while in use. 18.5" w x 11.5" d x 8" high. Shipping weight 6.0 lbs. (2.7 kg)

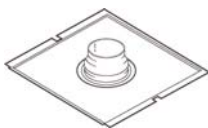
PURIFIER® CLASS I & HEPA FILTERED SAFETY ENCLOSURES

ACCESSORIES & DUCTWORK




Remote Blowers

Coated steel blowers with belt-drive motors. See page 28 for more information.



Canopy Connections

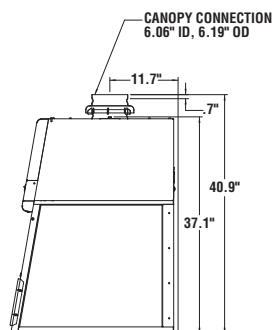Connection allows the Enclosure to be canopy-ducted to the outside for additional protection. The canopy's air gap balances the air by allowing room air to enter the ductwork and mix with the filtered air being exhausted from the Enclosure. Outlet accepts 6" nominal diameter ductwork. **Ductwork and house exhaust or Remote Blower are required (not included).**



Catalog #	Description	For use with:	Exhaust Volume	Shipping Weight
3924400	2-Ft Canopy Connection	2-Ft Purifier Class I & HEPA Filtered Enclosures	110 to 155 CFM	15 lbs. (6.8 kg)
3924401	3-Ft Canopy Connection	3-Ft Purifier Class I & HEPA Filtered Enclosures	165 to 230 CFM	18 lbs. (8.2 kg)
3924402	4-Ft Canopy Connection	4-Ft Purifier Class I & HEPA Filtered Enclosures	220 to 310 CFM	21 lbs. (9.5 kg)

Thermoplastic Duct

PVC exhaust duct in 10' lengths. A Female Duct Coupling and solvent cement are required to join two sections.

Catalog #	Nominal Dia.	Actual OD	Actual ID	Shipping Wt.
4708600	6"	6.625"	6.25"	25 lbs. (11.3 kg)
4718900	8"	8.625"	8.25"	35 lbs. (15.9 kg)

Female Duct Coupling

PVC coupling joins two sections of Thermoplastic Duct.

Catalog #	Nominal Dia.	Equivalent Resistance*	Shipping Wt.
4708900	6"	0	4 lbs. (1.8 kg)

Male Duct Coupling

PVC coupling joins two Duct Reducers and Duct Reducer to Blower.

Catalog #	Nominal Dia.	Equivalent Resistance*	Shipping Wt.
4719900	8"	0	4 lbs. (1.8 kg)
7027800	10"	0	5 lbs. (2.3 kg)

90° Elbow

PVC Elbow has belled end connections to receive Thermoplastic Duct or Male Duct Coupling directly.

Catalog #	Nominal Dia.	Equivalent Resistance*	Shipping Wt.
4708700	6"	12 ft.	8 lbs. (3.6 kg)

*Equivalent resistance is measured in feet of straight duct.

Flexible Duct Connection

Reduces vibration between the blower and PVC ductwork. Supplied with two clamps.

Catalog #	Nominal Dia.	Shipping Wt.
4861800	7" for use with 6" fittings	8 lbs. (3.6 kg)

Exhaust Dampers

Catalog #	Nominal Diameter	Material	Equivalent Resistance*	Shipping Weight
4724200	6"	PVC	1 ft.	1 lb. (0.5 kg)
3924000	6"	Powder-coated steel	1 ft.	1 lb. (0.5 kg)

Duct Reducers

PVC reducers connect duct and male couplings of different diameters.

Catalog #	Nominal Dia.	Equivalent Resistance*	Shipping Wt.
5605900	6" x 8"	0	2 lbs. (0.9 kg)
5606000	8" x 10"	0	5 lbs. (2.3 kg)

Coated-Coated Steel Blower Transition Adapter


Catalog #	Nominal Dia.	Equivalent Resistance*	Shipping Wt.
4722400	8"	1 ft.	3 lbs. (1.4 kg)

Zero Pressure Weathercap

PVC Weathercap fits atop standard PVC duct, permits vertical discharge of effluent air above roofline for dispersion away from building.

Catalog #	Nominal Diameter	Equivalent Resistance*	Height	Shipping Weight
4722300	8"	5 ft.	40"	25 lbs. (11.3 kg)

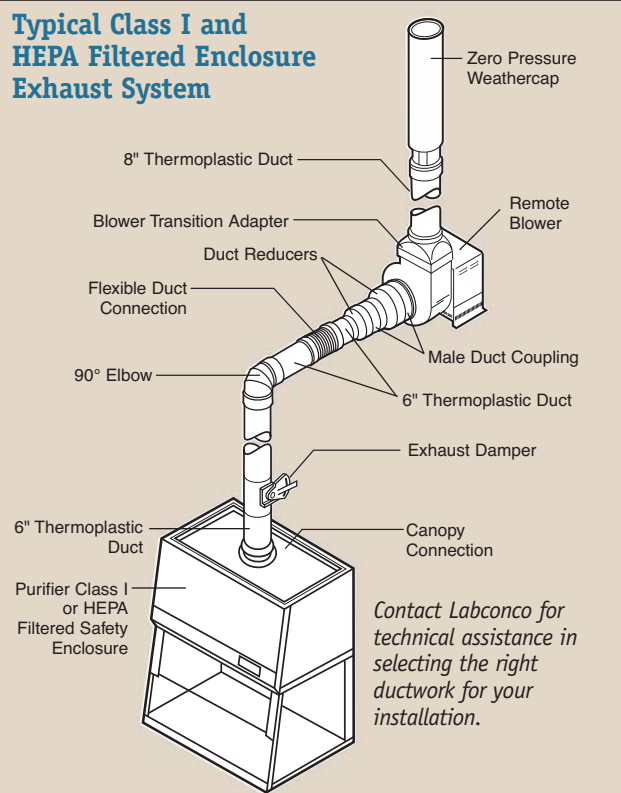Typical Class I and HEPA Filtered Enclosure Exhaust System



PURIFIER® HORIZONTAL CLEAN BENCHES

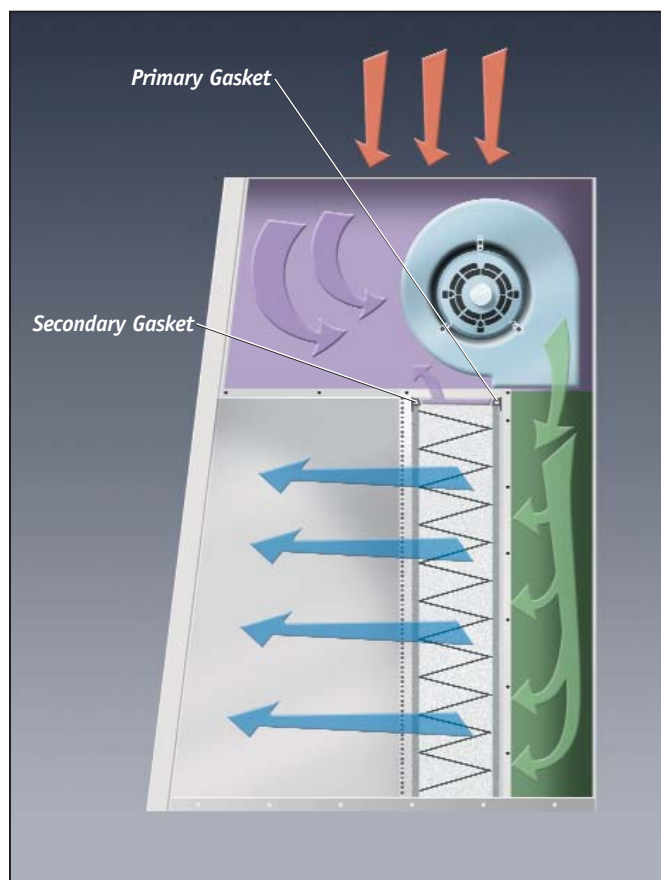
I N T R O D U C T I O N

The Purifier Horizontal Clean Bench protects your work. It is the ISO Class 5 work station of choice to protect materials that do not generate harmful aerosols or vapors, yet require a particulate-free environment. Suitable applications for the bench include plant tissue culture research, media preparation, electronic part inspection, syringe filling, medical device assembly and parenteral drug formulation.

Because air from the work area is dispersed directly into the laboratory, Purifier Horizontal Clean Benches should never be used in conjunction with biohazardous material, toxins or radionuclides. You and your safety officer must carefully assess the risk associated with any operation performed in a clean bench. Purifier Horizontal Clean Benches are offered in 3', 4' and 6' widths. Contact Labconco for information on additional widths.



Purifier Horizontal Clean Bench 3612504 on Telescoping Base Stand 3746702 with Ergonomic Chair 3744000.



Room Air

HEPA-Filtered Air

Unfiltered Air under Negative Pressure

Unfiltered Air under Positive Pressure

During operation, room air is drawn through a prefilter in the top of the bench to trap large particles. Then, the air flows through a 99.99% efficient HEPA filter and is projected horizontally across the entire work area. This HEPA-filtered air minimizes cross contamination in the work area and provides a particulate-free work environment.

Unique to the Purifier Horizontal Clean Bench is its double HEPA filter seal. The primary gasket seals the upstream side of the filter to prevent unfiltered air from leaking into the work area. The secondary gasket at the front of the HEPA filter seals the downstream side of the filter frame. The area between these two gaskets is maintained under negative pressure. Should a leak occur at the primary HEPA filter gasket, the contaminated air is immediately recaptured and refiltered.

PURIFIER® HORIZONTAL CLEAN BENCHES

FEATURES & BENEFITS

Replaceable prefilters trap large particles, extend HEPA filter life.

99.99% efficient HEPA filter is industry standard size for economical replacement.

Low vibration, variable speed motor/blower 1/3 hp on 3-foot models, 1/2 hp on 4-foot models, and two 1/4 hp motors on 6-foot models.

Minihelic II pressure gauge displays differential pressure across the HEPA filter.

Particle-free ISO Class 5 air. Particle counts in the work area conform to ISO 14644-1 and 2, providing ISO Class 5 conditions (formerly Class 100). This ensures fewer than 3520 particles 0.5 µm or larger per cubic meter of air.

Full one year warranty on parts and service.

Optional Vibration Isolation Table. See page 11 for ordering information.

Glare-free fluorescent lighting.

CE Mark. All 230 volt models conform to the following CE (European Community requirements for electrical safety and electromagnetic compatibility).

Optional UV Screen acts as a barrier when bench is not in use. See page 11 for ordering information.

Accessory Telescoping Base Stand. See pages 24 and 25 for ordering information.

ETL listed. All 115 volt, 60 Hz models carry the ETL mark signifying that they conform to UL Standard 3101-1/61010-1 in the U.S. and are certified to CAN/CSA C22.2 No. 1010.1 in Canada.



Accessory Hydraulic Lift Base Stand. See pages 26 and 27 for ordering information.




 Exclusive feature

Front-mounted light and blower switches.

 **Angled profile design.** Front panel is free of protrusions that could interfere with visibility or obstruct taller users.



Polarized and grounded electrical outlet(s).

 **Double-gasket, negative pressure leak protection.**

Optional service fixture available factory-installed or as a kit for user installation. See page 11 for kit information.

Dark gray, melamine-laminated hard board work surface.

Optional Type 304 brushed stainless steel work surface and sides. Contact Labconco for ordering information.



Optional factory-installed 254 nm UV light for secondary decontamination while the bench is not in use. UV light switch is safety interlocked with fluorescent light and blower switches to help prevent inadvertent UV exposure while the bench is in operation.

PURIFIER® HORIZONTAL CLEAN BENCHES

ORDERING INFORMATION

Standard Features

- 99.99% efficient HEPA filter
- Prefilter
- Class 5 conditions per ISO 14644-1 and 2 (formerly Class 100)
- Minihelic II pressure gauge
- Double-gasket, negative pressure HEPA filter leak protection
- Front-mounted light and blower switches
- Variable speed motor/blower(s) with vibration isolation pads
- Fluorescent lighting
- Polarized and grounded electrical outlet(s)
- Glacier white epoxy-coated steel frame and exterior panels
- Dark gray melamine-laminated hard board work surface
- Angled front profile
- ETL listing, UL 3101-1/61010-1 and CAN/CSA C22.2 No. 1010.1 conformity (115 volt models)
- CE Conformity Marking (230 volt models)
- One year warranty on parts and service

Standard Option Package

- Factory-installed, 254 nm UV lamp with interlocking safety switch allowing operation only when blower and fluorescent light are off
- Factory-installed, chrome-plated, forged brass service fixture(s) with quarter turn handle

Required Accessory

- Supporting base. Place on existing casework or see pages 24-27.



Purifier Horizontal Clean Bench 3610004 on Telescoping Base Stand 3746704.

See page 11 for dimensional drawings.

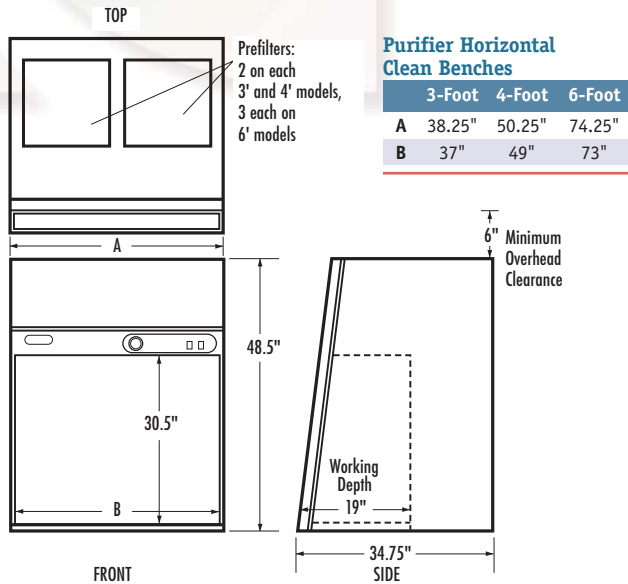
 Exclusive feature

Catalog #	Nominal Width**	Electrical Requirements	Power Cord & Plug	Light(s)	Electrical Outlets	Service Fixture(s)	Shipping Weight
3600000	3 feet	115 volts, 60 Hz, 12 amps	115 volts, 15 amps	Fluorescent	(1) 115 volt	—	305 lbs. (138 kg)
3600004	3 feet	115 volts, 60 Hz, 12 amps	115 volts, 15 amps	Fluorescent Ultraviolet	(1) 115 volt	1	305 lbs. (138 kg)
3600020*	3 feet	230 volts, 50 Hz, 7 amps	230 volts, 15 amps no plug	Fluorescent	(1) 230 volt (IEC 320)	—	305 lbs. (138 kg)
3600024*	3 feet	230 volts, 50 Hz, 7 amps	230 volts, 15 amps no plug	Fluorescent Ultraviolet	(1) 230 volt (IEC 320)	1	305 lbs. (138 kg)
3612500	4 feet	115 volts, 60 Hz, 12 amps	115 volts, 15 amps	Fluorescent	(1) 115 volt	—	370 lbs. (168 kg)
3612504	4 feet	115 volts, 60 Hz, 12 amps	115 volts, 15 amps	Fluorescent Ultraviolet	(1) 115 volt	1	370 lbs. (168 kg)
3612520*	4 feet	230 volts, 50 Hz, 7 amps	230 volts, 15 amps no plug	Fluorescent	(1) 230 volt (IEC 320)	—	370 lbs. (168 kg)
3612524*	4 feet	230 volts, 50 Hz, 7 amps	230 volts, 15 amps no plug	Fluorescent Ultraviolet	(1) 230 volt (IEC 320)	1	370 lbs. (168 kg)
3610000	6 feet	115 volts, 60 Hz, 16 amps	115 volts, 20 amps	Fluorescent	(2) 115 volt	—	515 lbs. (234 kg)
3610004	6 feet	115 volts, 60 Hz, 16 amps	115 volts, 20 amps	Fluorescent Ultraviolet	(2) 115 volt	2	515 lbs. (234 kg)
3610020*	6 feet	230 volts, 50 Hz, 8 amps	230 volts, 15 amps no plug	Fluorescent	(2) 230 volt (IEC 320)	—	515 lbs. (234 kg)
3610024*	6 feet	230 volts, 50 Hz, 8 amps	230 volts, 15 amps no plug	Fluorescent Ultraviolet	(2) 230 volt (IEC 320)	2	515 lbs. (234 kg)

*International electrical configuration **Contact Labconco for additional widths.

PURIFIER® HORIZONTAL CLEAN BENCHES

DIMENSIONS & ACCESSORIES



UV Screen Kits

Provide a barrier to ultraviolet light while UV light is on. Valance mounts above opening and hides retracted screen when not in use. Include screen, valance and mounting brackets. **Installation required.**

Catalog #	For use with:	Shipping Weight
3617000	3-Ft Purifier Clean Benches	10 lbs. (4.5 kg)
3617001	4-Ft Purifier Clean Benches	12 lbs. (5.4 kg)
3617002	6-Ft Purifier Clean Benches	14 lbs. (6.4 kg)



Replacement Prefilters

Labconco recommends replacing pre-filters every three months. One each. Shipping weight 1.0 lb. (0.4 kg)

Catalog #	For use with:	# Required
3768900	3-Ft Purifier Clean Benches	2
3768901	4-Ft Purifier Clean Benches	2
3768901	6-Ft Purifier Clean Benches	3



Telescoping Base Stands

Epoxy-coated steel frame. Adjust in 1" increments to provide a working height from 28.75 to 34.75". Available with fixed feet or casters. See pages 24 and 25 for more information.



Hydraulic Lift Base Stands

Epoxy-coated steel, with infinite height adjustment using push button or hand crank. See pages 26 and 27 for more information.



3697500 IV Bar Kit

Supports intravenous solution bottles and bags. Kit includes IV bar, mounting hardware, four hangers and instructions for installation. Three-foot and four-foot Purifier Clean Benches accommodate one IV bar. Six-foot benches accommodate two. Shipping weight 5.0 lbs. (2.3 kg)



3773700 Service Fixture Kit

Includes serrated hose tip valve with quarter turn control handle, hardware and instructions for plumbing to services. Up to two fixtures may be mounted on the right hand side of the Purifier Horizontal Clean Bench. Shipping weight 4.0 lbs. (1.8 kg)



3618000 Vibration Isolation Table

Provides an isolated work surface for conducting procedures with vibration-sensitive equipment such as microscopes and balances. The table never makes contact with the clean bench so vibration from the motor/blower is not transmitted to the table. Epoxy-coated steel frame, adjusts in 1" increments to provide a working height from 29.5 to 36". ADA-compliant. Four leveling feet. Work surface is laminated hard board. 30.4" w x 18" d. Overall dimensions: 32.5" x 25" d x 29.5" to 36" high. Shipping weight 110.0 lbs. (49.9 kg)



3744000 Ergonomic Chair with Armrests

Lab chair has 6-way articulating seat and back control for personalized adjustment. Pneumatic mechanism adjusts seat height from 18.25" to 25.75". Shipping weight 35.0 lbs. (15.9 kg). See page 66 for more information.



3746000 Adjustable Footrest

Elevates feet and permits angle repositioning while in use. 18.5" w x 11.5" d x 8" high. Shipping weight 6.0 lbs. (2.7 kg)

PURIFIER® VERTICAL CLEAN BENCHES

I N T R O D U C T I O N

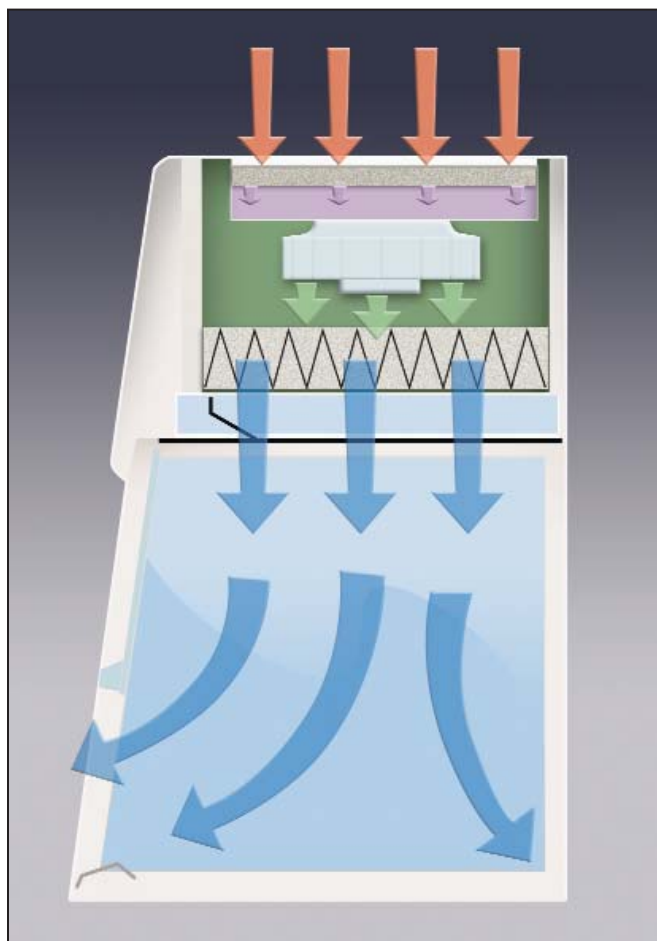
Purifier Vertical Clean Benches protect your work from particulate contamination. Using a 99.99% efficient supply HEPA filter, Purifier Vertical Clean Benches maintain ISO Class 5 conditions (formerly Class 100) in the work area. Appropriate applications for Clean Benches include plant tissue culture, media plate preparation, electronics inspection, medical device assembly and IV solution preparation.

Because Purifier Vertical Clean Benches do not provide protection to the user, they should not be used in conjunction with biohazardous material, toxins or radionuclides. You and your safety officer must carefully assess the risk associated with any operation performed in a clean bench.


Purifier Vertical Clean Benches are offered in 2', 3' and 4' widths. Built-in options include a UV light for secondary decontamination, bi-fold doors for complete closure and a Guardian Airflow Monitor to monitor downflow velocity and HEPA filter loading.



Purifier Vertical Clean Bench 3970200 on Solid Epoxy Dished Work Surface 3909900 and Telescoping Base Stand 3746710 with Ergonomic Chair 3744000 and Adjustable Footrest 3746000.



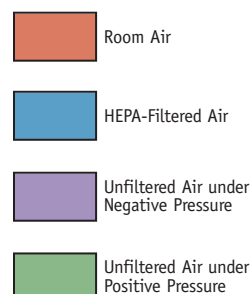During operation, room air is drawn through a prefilter in the top of the cabinet to trap large particles. Then, the air flows through a 99.99% efficient HEPA filter and is projected vertically over the work area. This HEPA-filtered air prevents contaminants from entering the work area and provides a particulate-free work environment that minimizes cross contamination.



PURIFIER® VERTICAL CLEAN BENCHES

FEATURES & BENEFITS

Particulate-free ISO Class 5 air. Particle counts in the work area conform to ISO 14644-1 and 2, providing ISO Class 5 conditions (formerly Class 100 per Federal Standard 209E), ensuring fewer than 3520 particles 0.5 µm or larger per cubic meter of air.

Front-mounted Minihelic pressure gauge indicates the total system pressure across the HEPA filter and pressurized plenum to help determine HEPA filter loading.

Disposable prefilter traps large particles to extend the life of the HEPA filter.

Optional 254 nm ultraviolet lamp provides secondary decontamination while the bench is not in use. The UV light does not penetrate the plane of the sash and air foil.

HEPA filter is rated to remove 99.99% of all particles 0.3 micron in size. It is industry standard size for economical replacement.

Full one year warranty on parts and service.

Air diffuser protects the HEPA filter from possible damage from direct contact with the operator and apparatus inside the enclosure. It distributes air evenly across the work area to optimize airflow.

Variable speed blower with solid state control may be accessed by a certified technician to validate and adjust downflow velocity.

Light and blower switches are located within easy reach of the operator. On models with UV light, a single 3-way switch powers the fluorescent and UV light, but only one at a time, helping to protect the operator from inadvertent UV exposure.

Tempered safety glass sash is sloped 10° for less glare and closer, more comfortable viewing than vertical sashes offer. The glass is UV opaque and resistant to UV irradiation. Wiping seals ensure ISO Class 5 conditions inside the enclosure. The sash has a spring-loaded latch so that it may be pivoted upward and locked open for loading and cleaning.

CE mark. All 230 volt, 50 Hz models conform to the CE (European Community) requirements for electrical safety and electromagnetic compatibility.

ETL listed. All 115 volt, 60 Hz models carry the ETL mark signifying that they are certified to UL Standard 3101-1/61010-1 and CAN/CSA C22.2 No. 1010.1.

Two convenient utility ports with iris openings allow passage of tubing and electrical cords from equipment inside the enclosure through the back for connection to services. Plugs are provided to seal the ports when not in use.

Optional bi-fold doors, found on models with UV lamps, allow the sash opening to be closed while UV light is on and prevent contaminants from entering when the clean bench is not in use.

Accessory Solid Epoxy Dished Work Surface. See page 55 for ordering information.

Glare-free fluorescent lighting. The lamp is located above the work area, out of contact with the filtered air.

Optional built-in Guardian™ Airflow Monitor alerts the user to downflow velocities outside of preset range and HEPA filter conditions.



Exclusive feature



PURIFIER® VERTICAL CLEAN BENCHES

ORDERING INFORMATION

Standard Features

- Nominal downflow velocity of 40-55 fpm
- Supply HEPA filter, 99.99% efficient on particles 0.3 micron
- ISO Class 5 air (formerly Class 100 per Federal Standard 209E)
- Disposable prefilter
- Variable speed built-in blower with motorized impeller and solid state control rated for 5 amps
- Minihelic pressure gauge
- Air diffuser
- UV-resistant, glacier white and gray epoxy-coated aluminum and steel construction
- UV-opaque and resistant 1/4" thick tempered safety glass sash and sides
- Fluorescent lamp
- Light and blower switches
- 9.4" high sash opening
- Ergonomic 10° angled and hinged sash that pivots upward, locks to a loading height of 20" and has wiping seals to promote ISO Class 5 conditions
- Two utility ports with iris openings and plugs
- ETL listing, UL 3101-1/61010-1 and CAN/CSA C222.2 No. 1010.1 conformity (115 volt models)
- CE Conformity Marking (230 volt models)
- One year warranty on parts and service

Standard Options

- Guardian Airflow Monitor that continuously monitors downflow airflow and has two LED lights to indicate when airflow is within or outside set point range. Set point is factory calibrated for 25-35 fpm downflow velocity. The "Safe" green LED glows when airflow is within set point range. The "Alert" red LED glows and an audible alarm sounds when airflow is outside set point range. The audible alarm may be muted by pressing the "Silence Alarm" switch.
- 254 nm UV lamp with 3-way switch (fluorescent, UV and off) and bi-fold doors that close the sash opening.

Required Accessories

- Supporting work surface. Place on existing casework or see page 19.
- Supporting base. Place on existing casework or see pages 24-27.

See page 19 for dimensional drawings.

■ Exclusive feature



Purifier Vertical Clean Bench 3970200 on Solid Epoxy Dished Work Surface 3909900 and Telescoping Base Stand 3746710.

PURIFIER® VERTICAL CLEAN BENCHES

ORDERING INFORMATION

Catalog #	Nominal Width	Electrical Requirements	UV Lamp/ Bi-Fold Doors	Airflow Monitor	Power Cord & Plug	Shipping Weight
3970200	2 feet	115 volts, 60 Hz, 3 amps			115 volts, 15 amps	125 lbs. (57 kg)
3970201	2 feet	115 volts, 60 Hz, 3 amps	•		115 volts, 15 amps	125 lbs. (57 kg)
3970203	2 feet	115 volts, 60 Hz, 3 amps		•	115 volts, 15 amps	125 lbs. (57 kg)
3970204	2 feet	115 volts, 60 Hz, 3 amps	•	•	115 volts, 15 amps	125 lbs. (57 kg)
3970220*	2 feet	230 volts, 50 Hz, 2 amps			230 volts, 10 amps, no plug	125 lbs. (57 kg)
3970221*	2 feet	230 volts, 50 Hz, 2 amps	•		230 volts, 10 amps, no plug	125 lbs. (57 kg)
3970223*	2 feet	230 volts, 50 Hz, 2 amps		•	230 volts, 10 amps, no plug	125 lbs. (57 kg)
3970224*	2 feet	230 volts, 50 Hz, 2 amps	•	•	230 volts, 10 amps, no plug	125 lbs. (57 kg)
3970300	3 feet	115 volts, 60 Hz, 5 amps			115 volts, 15 amps	160 lbs. (73 kg)
3970301	3 feet	115 volts, 60 Hz, 5 amps	•		115 volts, 15 amps	160 lbs. (73 kg)
3970303	3 feet	115 volts, 60 Hz, 5 amps		•	115 volts, 15 amps	160 lbs. (73 kg)
3970304	3 feet	115 volts, 60 Hz, 5 amps	•	•	115 volts, 15 amps	160 lbs. (73 kg)
3970320*	3 feet	230 volts, 50 Hz, 3 amps			230 volts, 10 amps, no plug	160 lbs. (73 kg)
3970321*	3 feet	230 volts, 50 Hz, 3 amps	•		230 volts, 10 amps, no plug	160 lbs. (73 kg)
3970323*	3 feet	230 volts, 50 Hz, 3 amps		•	230 volts, 10 amps, no plug	160 lbs. (73 kg)
3970324*	3 feet	230 volts, 50 Hz, 3 amps	•	•	230 volts, 10 amps, no plug	160 lbs. (73 kg)
3970400	4 feet	115 volts, 60 Hz, 5 amps			115 volts, 15 amps	195 lbs. (88 kg)
3970401	4 feet	115 volts, 60 Hz, 5 amps	•		115 volts, 15 amps	195 lbs. (88 kg)
3970403	4 feet	115 volts, 60 Hz, 5 amps		•	115 volts, 15 amps	195 lbs. (88 kg)
3970404	4 feet	115 volts, 60 Hz, 5 amps	•	•	115 volts, 15 amps	195 lbs. (88 kg)
3970420*	4 feet	230 volts, 50 Hz, 3 amps			230 volts, 10 amps, no plug	195 lbs. (88 kg)
3970421*	4 feet	230 volts, 50 Hz, 3 amps	•		230 volts, 10 amps, no plug	195 lbs. (88 kg)
3970423*	4 feet	230 volts, 50 Hz, 3 amps		•	230 volts, 10 amps, no plug	195 lbs. (88 kg)
3970424*	4 feet	230 volts, 50 Hz, 3 amps	•	•	230 volts, 10 amps, no plug	195 lbs. (88 kg)

*International electrical configuration

PURIFIER® FILTERED PCR ENCLOSURES

I N T R O D U C T I O N

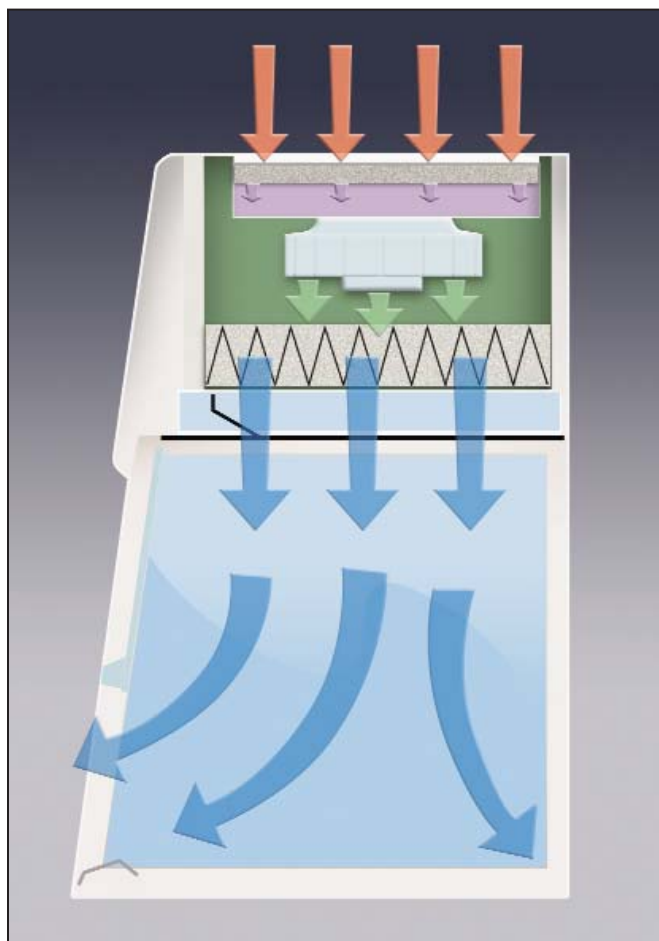
Purifier Filtered PCR Enclosures are vertical clean benches that protect your work. Available in 2', 3' and 4' widths, they provide a controlled environment in which to perform polymerase chain reaction experiments. ISO Class 5 air is constantly projected vertically down over the work area, minimizing the risk of cross contamination of samples.

Purifier Filtered PCR Enclosures do not provide protection to the user and should not be used in conjunction with biohazardous material, toxins or radionuclides. You and your safety officer must carefully assess the risks associated with any operation performed in the PCR enclosure.


The variable digital timer provides UV light exposure continuously or in user-selected increments of 5, 10, 15, 30, 60, 120 or 240 minutes to deactivate DNA and RNA contaminants. When the timer expires, the UV light automatically switches off in preparation for the next experiment.



Purifier Filtered PCR Enclosure 3970305 on Solid Epoxy Dished Work Surface 3909901 and Telescoping Base Stand 3746711 with Ergonomic Chair 3744000 and Adjustable Footrest 3746000.



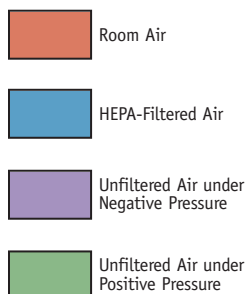During operation, room air is drawn through a prefilter in the top of the cabinet to trap large particles. Then, the air flows through a 99.99% efficient HEPA filter and is projected vertically over the work area. This HEPA-filtered air prevents contaminants from entering the work area and provides a particulate-free work environment that minimizes cross contamination.



PURIFIER® FILTERED PCR ENCLOSURES

FEATURES & BENEFITS

Particulate-free ISO Class 5 air. Particle counts in the work area conform to ISO 14644-1 and 2, providing ISO Class 5 conditions (formerly Class 100 per Federal Standard 209E), ensuring fewer than 3520 particles 0.5 µm or larger per cubic meter of air.

Front-mounted Minihelic pressure gauge indicates the total system pressure across the HEPA filter and pressurized plenum to help determine HEPA filter loading.

ETL listed. All 115 volt, 60 Hz models carry the ETL mark signifying that they are certified to UL Standard 3101-1/61010-1 and CAN/CSA C22.2 No. 1010.1.

CE mark. All 230 volt, 50 Hz models conform to the CE (European Community) requirements for electrical safety and electromagnetic compatibility.

Two convenient utility ports with iris openings allow passage of tubing and electrical cords from equipment inside the enclosure through the back for connection to services. Plugs are provided to seal the ports when not in use.

Bi-fold doors allow the sash opening to be closed while the UV light is on and prevent contaminants from entering when the enclosure is not in use.

Variable digital timer provides timed UV light exposure. From the front-mounted control panel, the operator may select from eight settings in minutes of 5, 10, 15, 30, 60, 120, 240 and continuous. When the UV light switch is activated, the previously selected time setting is automatically initiated or the user may select a different setting. An amber LED indicator blinks for the duration of the setting.

Disposable prefilter traps large particles to extend the life of the HEPA filter.

The 254 nm ultraviolet lamp denatures stray DNA strands and provides secondary decontamination while the enclosure is not in use.

HEPA filter is rated to remove 99.99% of all particles 0.3 micron in size. It is industry standard size for economical replacement.

Full one year warranty on parts and service.

Air diffuser protects the HEPA filter from possible damage from direct contact with the operator and apparatus inside the enclosure. It distributes air evenly across the work area to optimize airflow.

Variable speed blower with solid state control may be accessed by a certified technician to validate and adjust downflow velocity.

Glare-free fluorescent lighting. The lamp is located above the work area, out of contact with the filtered air.

Light and blower switches are located within easy reach of the operator. The 3-way light switch powers the fluorescent and UV light, but only one at a time, helping to protect the operator from inadvertent UV exposure.

Tempered safety glass sash is sloped 10° for less glare and closer, more comfortable viewing than vertical sashes offer. The glass is UV opaque and resistant to UV irradiation. Wiping seals ensure ISO Class 5 conditions inside the enclosure. The sash has a spring-loaded latch so that it may be pivoted upward and locked open for loading and cleaning.



Accessory Solid Epoxy Dished Work Surface. See page 55 for ordering information.



Optional built-in Guardian™ Airflow Monitor alerts the user to downflow velocities outside of preset range and HEPA filter conditions.



PURIFIER® FILTERED PCR ENCLOSURES

ORDERING INFORMATION

Standard Features

- Nominal downflow velocity of 40-55 fpm
- Supply HEPA filter, 99.99% efficient on particles 0.3 micron
- ISO Class 5 air (formerly Class 100 per Federal Standard 209E)
- Disposable prefilter
- Variable speed built-in blower with motorized impeller and solid state control rated for 5 amps
- Minihelic pressure gauge
- Air diffuser
- UV-resistant, glacier white and gray epoxy-coated aluminum and steel construction
- UV-opaque and resistant 1/4" thick tempered safety glass sash and sides
- 9.4" high sash opening
- Fluorescent lamp
- 254 nm UV lamp
- Variable digital timer for UV exposure with eight settings in minutes of 5, 10, 15, 30, 60, 120, 240 and continuous
- Blower and 3-way light (fluorescent, UV and off) switches
- Ergonomic 10° angled and hinged sash that pivots upward, locks to a loading height of 20" and has wiping seals to promote ISO Class 5 conditions
- Bi-fold doors that close the sash opening
- Two utility ports with iris openings and plugs
- ETL listing, UL 3101-1/61010-1 and CAN/CSA C222.2 No. 1010.1 conformity (115 volt models)
- CE Conformity Marking (230 volt models)
- One year warranty on parts and service



Purifier Filtered PCR Enclosure 3970305 on Solid Epoxy Dished Work Surface 3909901 and Telescoping Base Stand 3746711.

Standard Option

- Guardian Airflow Monitor that continuously monitors down-flow airflow and has two LED lights to indicate when airflow is within or outside set point range. Set point is factory calibrated for 25-35 fpm downflow velocity. The "Safe" green LED glows when airflow is within set point range. The "Alert" red LED glows and an audible alarm sounds when airflow is outside set point range. The audible alarm may be muted by pressing the "Silence Alarm" switch.

Required Accessories

- Supporting work surface. Place on existing casework or see page 19.
- Supporting base. Place on existing casework or see pages 24-27.

See page 19 for dimensional drawings.


Catalog #	Nominal Width	Electrical Requirements	Variable Timed UV Lamp/Bi-Fold Doors	Airflow Monitor	Power Cord & Plug	Shipping Weight
3970202	2 feet	115 volts, 60 Hz, 3 amps	•		115 volts, 15 amps	125 lbs. (57 kg)
3970205	2 feet	115 volts, 60 Hz, 3 amps	•	•	115 volts, 15 amps	125 lbs. (57 kg)
3970222*	2 feet	230 volts, 50 Hz, 2 amps	•		230 volts, 10 amps, no plug	125 lbs. (57 kg)
3970225*	2 feet	230 volts, 50 Hz, 2 amps	•	•	230 volts, 10 amps, no plug	125 lbs. (57 kg)
3970302	3 feet	115 volts, 60 Hz, 3 amps	•		115 volts, 15 amps	160 lbs. (73 kg)
3970305	3 feet	115 volts, 60 Hz, 3 amps	•	•	115 volts, 15 amps	160 lbs. (73 kg)
3970322*	3 feet	230 volts, 50 Hz, 2 amps	•		230 volts, 10 amps, no plug	160 lbs. (73 kg)
3970325*	3 feet	230 volts, 50 Hz, 2 amps	•	•	230 volts, 10 amps, no plug	160 lbs. (73 kg)
3970402	4 feet	115 volts, 60 Hz, 3 amps	•		115 volts, 15 amps	195 lbs. (88 kg)
3970405	4 feet	115 volts, 60 Hz, 3 amps	•	•	115 volts, 15 amps	195 lbs. (88 kg)
3970422*	4 feet	230 volts, 50 Hz, 2 amps	•		230 volts, 10 amps, no plug	195 lbs. (88 kg)
3970425*	4 feet	230 volts, 50 Hz, 2 amps	•	•	230 volts, 10 amps, no plug	195 lbs. (88 kg)

■ Exclusive feature


* International electrical configuration

PURIFIER® VERTICAL CLEAN BENCHES & FILTERED PCR ENCLOSURES

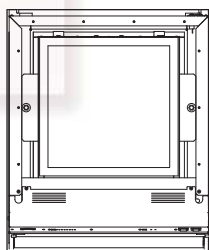
DIMENSIONS & ACCESSORIES



TOP

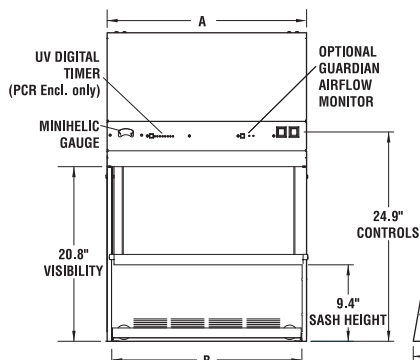

FRONT

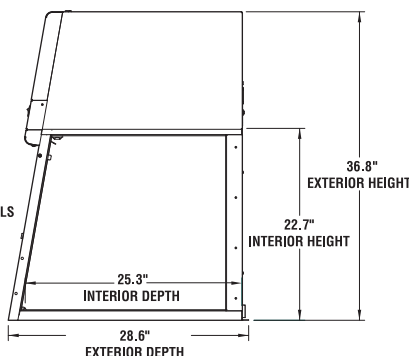

SIDE

Purifier Vertical Clean Benches & Filtered PCR Enclosures

	2-Foot	3-Foot	4-Foot
A	24.0"	36.0"	48.0"
B	22.8"	34.8"	46.8"



Replacement Disposable Prefilters

Corrugated paper prefilters trap large particles to extend HEPA filter life. Labconco recommends replacing the prefilter every three months. One each.

Catalog #	For use with	Shipping Weight
3935700	2-Ft Vertical Clean Bench	1.0 lb. (0.5 kg)
3935701	3-Ft Vertical Clean Bench	1.0 lb. (0.5 kg)
3935702	4-Ft Vertical Clean Bench	1.0 lb. (0.5 kg)



IV Bars

Support intravenous solution bottles and bags. Include IV bar, mounting hardware, four hangers and installation instructions.

Catalog #	For use with	Shipping Weight
3721000	2-Ft Vertical Clean Bench	4.0 lbs. (1.8 kg)
3721001	3-Ft Vertical Clean Bench	6.0 lbs. (2.7 kg)
3721002	4-Ft Vertical Clean Bench	8.0 lbs. (3.6 kg)



Base Stands

Epoxy-coated steel. Telescoping Base Stands adjust in 1" increments to provide a working height from 30" to 36" and are available with fixed feet or casters. Electric and Manual Hydraulic Lift Base Stands have

infinite height adjustment using a push button or hand crank. See pages 24-27 for more information.



Solid Epoxy Dished Work Surfaces

Black chemical-resistant work surface is contoured to fit the dimensions of the Purifier Vertical Clean Benches and Filtered PCR Enclosures and to contain spills.

Catalog #	For use with	Dimensions	Shipping Weight
3909900	2-Ft Purifier Enclosure	24.0" w x 29.0" d x 1.0" h (61 cm x 73.7 cm x 2.5 cm)	50.0 lbs. (23 kg)
3909901	3-Ft Purifier Enclosure	36.0" w x 29.0" d x 1.0" h (91.4 cm x 73.7 cm x 2.5 cm)	75.0 lbs. (34 kg)
3909902	4-Ft Purifier Enclosure	48.0" w x 29.0" d x 1.0" h (122 cm x 73.7 cm x 2.5 cm)	100.0 lbs. (45 kg)



3744000 Ergonomic Chair with Armrests

Chair has 6-way articulating seat and back control for personalized adjustment. Pneumatic mechanism adjusts seat height from 18.25" to 25.75". Shipping weight 35.0 lbs. (15.9 kg). See page 28 for more information.



3746000 Adjustable Footrest

Elevates feet and permits angle repositioning while in use. 18.5" w x 11.5" d x 8" high. Shipping weight 6.0 lbs. (2.7 kg)

PURIFIER® NON-VENTILATED PCR & TISSUE CULTURE ENCLOSURES

FEATURES & BENEFITS

Purifier Non-Ventilated Enclosures provide a circulation-free work area designed to reduce cross contamination. These enclosures have an ultraviolet light that decontaminates surfaces and denatures trace DNA during periods of non-use. Bi-fold doors close the sash opening while the UV light is in use to prevent inadvertent personnel exposure and contaminants from entering the enclosure.

The Purifier Non-Ventilated PCR Enclosures, designed for conducting polymerase chain reaction procedures,

include a UV light timer that allows the user to select from eight time settings. The Purifier Tissue Culture Enclosures may be used for culture manipulation of non-hazardous samples. Because these enclosures do not provide protection to the user, they should not be used in conjunction with biohazardous material, toxins or radionuclides. You and your safety officer must carefully assess the risk associated with any operation performed in a Purifier Non-Ventilated Enclosure.

Full one year warranty on parts and service.

Glare-free fluorescent lighting provides illumination inside the enclosure.

The 254 nm ultraviolet lamp eliminates stray DNA and provides secondary decontamination while the enclosure is not in use.

Electrical receptacle located inside the enclosure is powered by the front-mounted auxiliary switch.

Light and auxiliary electrical receptacle switches are located within easy reach of the operator. The 3-way light switch powers the fluorescent and UV light, but only one at a time, helping to protect the operator from inadvertent UV exposure.



Tempered safety glass sash is sloped 20° for less glare and closer, more comfortable viewing than vertical sashes offer. The glass is UV opaque and resistant to UV irradiation. Wiping seals reduce cross contamination inside the enclosure. The sash has a spring-loaded latch so that it may be pivoted upward and locked open for loading and cleaning.

CE mark. All 230 volt, 50 Hz models conform to the CE (European Community) requirements for electrical safety and electromagnetic compatibility.

Two convenient utility ports with iris openings allow passage of tubing and electrical cords from equipment inside the enclosure through the back for connection to services. Plugs are provided to seal the ports when not in use.

Bi-fold doors allow the sash opening to be closed while the UV light is on and prevent contaminants from entering when the enclosure is not in use.

Accessory Solid Epoxy Dished Work Surface. See page 59 for ordering information.

ETL listed. All 115 volt, 60 Hz models carry the ETL mark signifying that they are certified to UL Standard 3101-1/61010-1 and CAN/CSA C22.2 No. 1010.1.

Variable digital timer, found on Purifier Non-Ventilated PCR Enclosures, provides timed UV light exposure. From the front-mounted control panel, the operator may select from eight settings in minutes of 5, 10, 15, 30, 60, 120, 240 and continuous. When the UV light switch is activated, the previously selected time setting is automatically initiated or the user may select a different setting. An amber LED indicator blinks for the duration of the setting.








 Exclusive feature

PURIFIER® NON-VENTILATED PCR ENCLOSURES

ORDERING INFORMATION

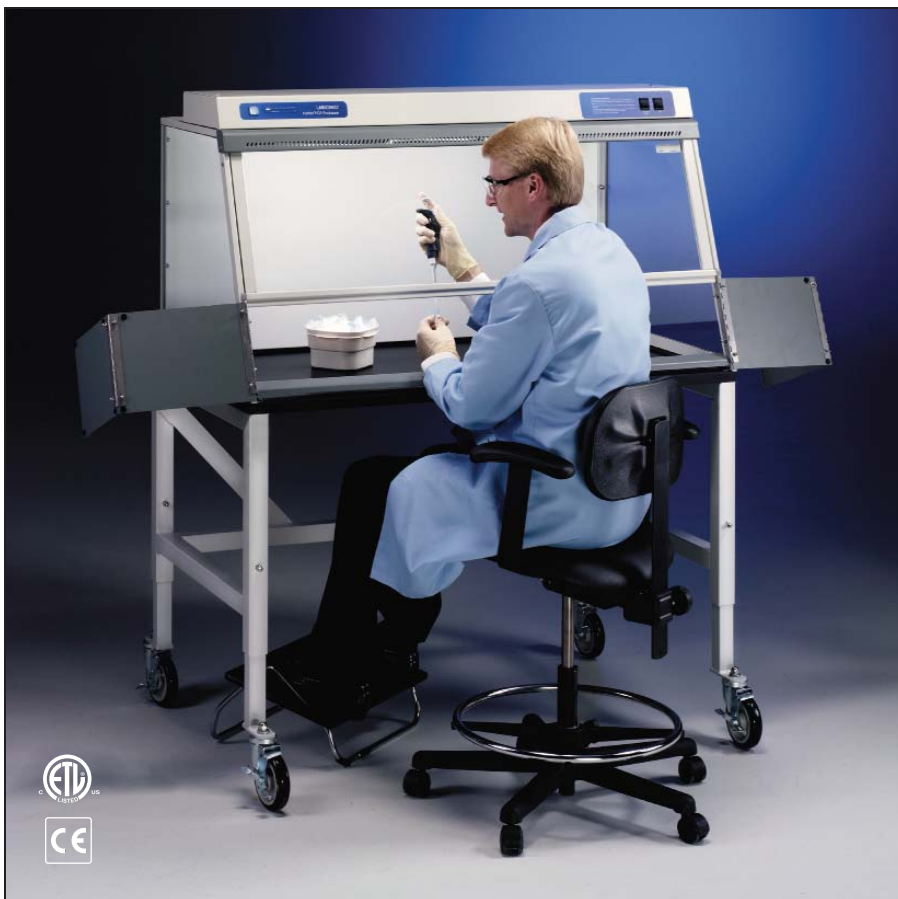
Standard Features

-  UV-resistant, glacier white and gray epoxy-coated aluminum and steel construction
-  UV-opaque and resistant, 1/4" thick tempered safety glass sash and sides
 - 7.8" high sash opening
-  Ergonomic 20° angled and hinged sash that pivots upward, locks to a loading height of 20" and has wiping seals to reduce cross contamination
-  Bi-fold doors that close the sash opening
 - Fluorescent lamp
 - 254 nm UV lamp
 - Interior electrical receptacle
-  Variable digital timer for UV exposure with eight settings in minutes of 5, 10, 15, 30, 60, 120, 240 and continuous
 - Front-mounted switches for 3-way light (fluorescent, UV and off) and auxiliary electrical receptacle
 - Two utility ports with iris openings and plugs
 - ETL listing, UL 3101-1/61010-1 and CAN/CSA C222.2 No. 1010.1 conformity (115 volt models)
 - CE Conformity Marking (230 volt models)
 - One year warranty on parts and service

Required Accessories

- Supporting work surface. Place on existing casework or see page 23.
- Supporting base. Place on existing casework or see pages 24 and 25.

See page 23 for dimensional drawings.



Purifier Non-Ventilated PCR Enclosure 3952400 on Solid Epoxy Dished Work Surface 3908402 and Telescoping Base Stand 3747012 with Ergonomic Chair 3744000 and Adjustable Footrest 3746000.





Catalog #	Nominal Width	Electrical Requirements	Variable Timed UV Lamp/Bi-Fold Doors	Power Cord & Plug	Shipping Weight
3952400	4 feet	115 volts, 60 Hz, 10 amps	•	115 volts, 15 amps	100 lbs. (45 kg)
3952420*	4 feet	230 volts, 50 Hz, 5 amps	•	230 volts, 10 amps, no plug	100 lbs. (45 kg)

*International electrical configuration

PURIFIER® NON-VENTILATED TISSUE CULTURE ENCLOSURES

ORDERING INFORMATION

Standard Features

-  UV-resistant, glacier white and gray epoxy-coated aluminum and steel construction
-  UV-opaque and resistant, 1/4" thick tempered safety glass sash and sides
 - 7.8" high sash opening
-  Ergonomic 20° angled and hinged sash that pivots upward, locks to a loading height of 20" and has wiping seals to reduce cross contamination
 - Fluorescent lamp
 - 254 nm UV lamp
 - Interior electrical receptacle
 - Front-mounted switches for 3-way light (fluorescent, UV and off) and auxiliary electrical receptacle
-  Bi-fold doors that close the sash opening
 - Two utility ports with iris openings and plugs
 - ETL listing, UL 3101-1/61010-1 and CAN/CSA C222.2 No. 1010.1 conformity (115 volt models)
 - CE Conformity Marking (230 volt models)
 - One year warranty on parts and service

Required Accessories

- Supporting work surface. Place on existing casework or see page 23.
- Supporting base. Place on existing casework or see pages 24 and 25.

See page 23 for dimensional drawings.




Purifier Tissue Culture Enclosure 3951400 on Solid Epoxy Dished Work Surface 3908402 and Telescoping Base Stand 3747012 with Ergonomic Chair 3744000 and Adjustable Footrest 3746000.

Catalog #	Nominal Width	Electrical Requirements	UV Lamp/ Bi-Fold Doors	Power Cord & Plug	Shipping Weight
3951400	4 feet	115 volts, 60 Hz, 10 amps	•	115 volts, 15 amps	100 lbs. (45 kg)
3951420*	4 feet	230 volts, 50 Hz, 5 amps	•	230 volts, 10 amps, no plug	100 lbs. (45 kg)

*International electrical configuration

PURIFIER® NON-VENTILATED PCR & TISSUE CULTURE ENCLOSURES

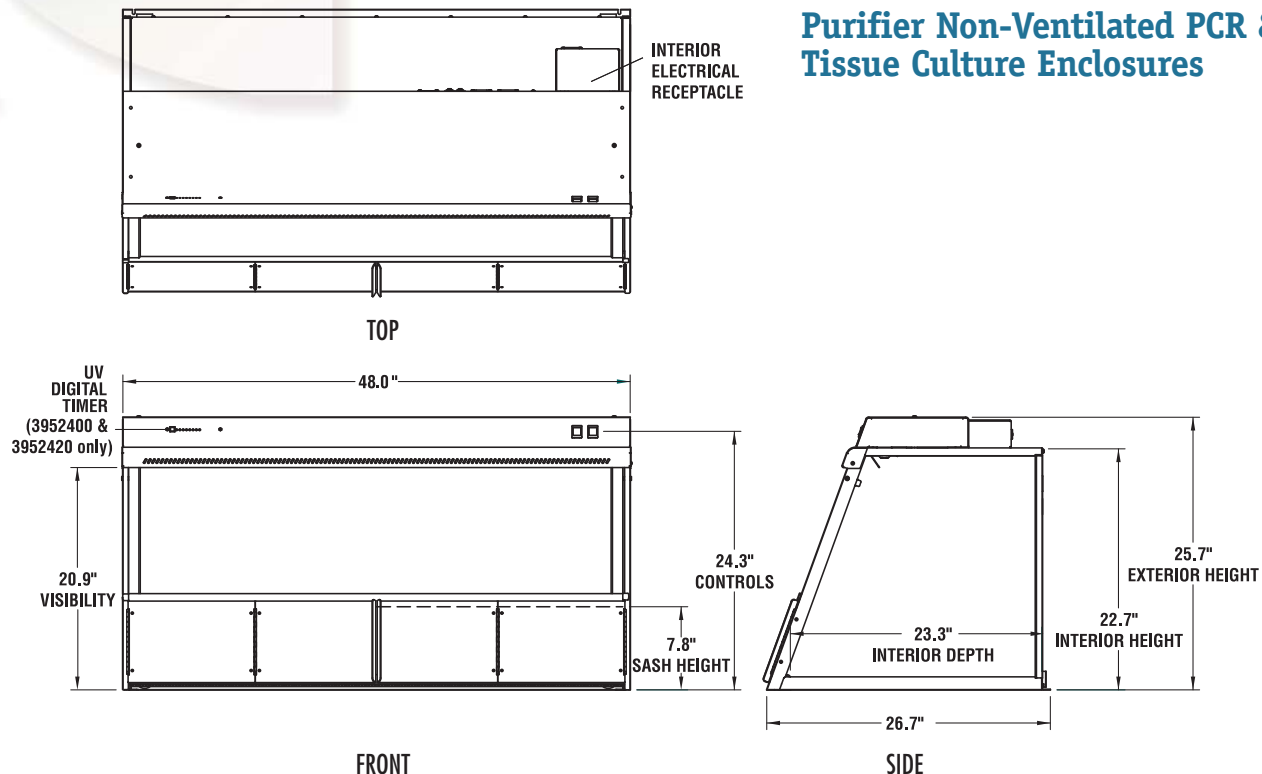D I M E N S I O N S & A C C E S S O R I E S



Solid Epoxy Dished Work Surface

Black chemical-resistant work surface is contoured to fit the dimensions of the Purifier Non-Ventilated PCR Enclosure and Tissue Culture Enclosure and to contain spills.

Catalog #	For use with	Dimensions	Shipping Weight
3908402	Non-Ventilated PCR & Tissue Culture Enclosures	48.0" w x 26.7" d x 1.0" h (122 cm x 67.8 cm x 2.5 cm)	96.0 lbs. (43.5 kg)



3744000 Ergonomic Chair with Armrests

Chair has 6-way articulating seat and back control for personalized adjustment. Pneumatic mechanism adjusts seat height from 18.25" to 25.75". Shipping weight 35.0 lbs. (15.9 kg). See page 28 for more information.



3746000 Adjustable Footrest

Elevates feet and permits angle repositioning while in use. 18.5" w x 11.5" d x 8" high. Shipping weight 6.0 lbs. (2.7 kg)



Telescoping Base Stands

Epoxy-coated steel. Adjust in 1" increments from 27.5" to 33.5" high. Available with fixed leveling feet or casters. See pages 24 and 25 for more information.

TELESCOPING BASE STANDS

These stands have telescoping legs so that they may be adjusted to seven height positions.* The stands are available with fixed feet or casters for mobility.

Standard Features

- Support any Purifier Delta Series Safety Cabinet, Purifier Clean Bench or Enclosure.
- Support loads up to 1000 pounds.
- ADA-compliant.
- Durable 1.75" tubular, epoxy-coated steel frame. Models 3' wide and larger have rear crossbar supports.
- Telescoping legs are adjustable to seven height positions in 1" increments from 27.5" to 33.5".*
- Available with four fixed leveling feet or 5" diameter, toe-locking, non-marking polyurethane casters with bearings for quiet operation.

*Height selection and adjustment are completed **before** the enclosure is installed atop the stand.

3730500 Caster Kit

For converting Telescoping Base Stands with Fixed Feet manufactured after 10/02. Kit includes four 5" diameter, non-marking polyurethane casters with toe locks. Load rating for each caster is 250 pounds. Casters have bearings for quiet operation. Casters replace the fixed feet included with Telescoping Base Stands. **Installation is required.** Shipping weight 10 lbs. (4.5 kg)




Telescoping Base Stand with Fixed Feet 3747010.

Inset: Telescoping Base Stand with Casters 3746701.




Catalog Numbers	Nominal Width	A	B	C
3746700, 3746710	2-Ft	24.0"	29.0"	8.2"
3746701, 3746711	3-Ft	36.0"	29.0"	8.2"
3747002, 3747012	4-Ft	48.0"	25.5"	8.2"
3746702, 3746712	4-Ft	48.0"	29.0"	8.2"
3746704, 3746714	6-Ft	72.0"	29.0"	8.2"

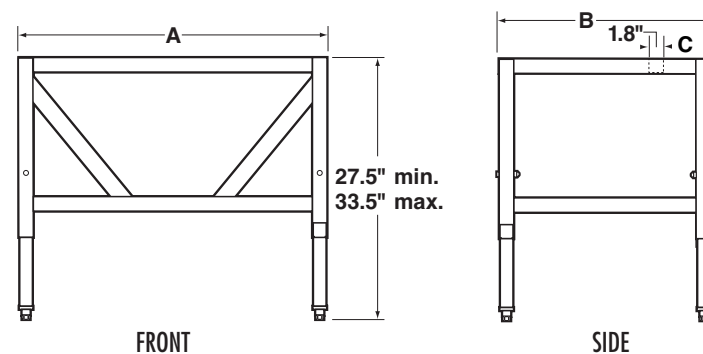With Fixed Feet



With Casters

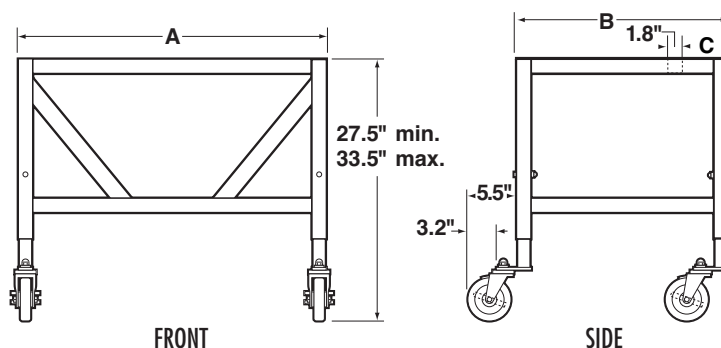

TELESCOPING BASE STANDS

Telescoping Base Stands with Fixed Feet

Nominal Width	Catalog #	Overall Width x Depth	Shipping Weight	For Use With:
2-Ft	3746700	24.0" w x 29.0" d	100 lbs. (45 kg)	Purifier Class I Enclosure, Purifier HEPA Filtered Enclosure, Purifier Vertical Clean Bench, Purifier Filtered PCR Enclosure
3-Ft	3746701	36.0" w x 29.0" d	100 lbs. (45 kg)	Purifier Class I Enclosure, Purifier HEPA Filtered Enclosure, Purifier Horizontal Clean Bench, Purifier Vertical Clean Bench, Purifier Filtered PCR Enclosure
4-Ft	3747002	48.0" w x 25.5" d	110 lbs. (50 kg)	Purifier Non-Ventilated PCR Enclosure, Purifier Tissue Culture Enclosure
4-Ft	3746702	48.0" w x 29.0" d	115 lbs. (52 kg)	Purifier Class I Enclosure, Purifier HEPA Filtered Enclosure, Purifier Horizontal Clean Bench, Purifier Vertical Clean Bench, Purifier Filtered PCR Enclosure
6-Ft	3746704	72.0" w x 29.0" d	125 lbs. (57 kg)	Purifier Horizontal Clean Bench

Telescoping Base Stands with 5" diameter Toe-Locking Casters*

Nominal Width	Catalog #	Overall Width x Depth	Shipping Weight	For Use With:
2-Ft	3746710	24.0" w x 29.0" d	100 lbs. (45 kg)	Purifier Class I Enclosure, Purifier HEPA Filtered Enclosure, Purifier Vertical Clean Bench, Purifier Filtered PCR Enclosure
3-Ft	3746711	36.0" w x 29.0" d	100 lbs. (45 kg)	Purifier Class I Enclosure, Purifier HEPA Filtered Enclosure, Purifier Horizontal Clean Bench, Purifier Vertical Clean Bench, Purifier Filtered PCR Enclosure
4-Ft	3747012	48.0" w x 25.5" d	110 lbs. (50 kg)	Purifier Non-Ventilated PCR Enclosure, Purifier Tissue Culture Enclosure
4-Ft	3746712	48.0" w x 29.0" d	115 lbs. (52 kg)	Purifier Class I Enclosure, Purifier HEPA Filtered Enclosure, Purifier Horizontal Clean Bench, Purifier Vertical Clean Bench, Purifier Filtered PCR Enclosure
6-Ft	3746714	72.0" w x 29.0" d	125 lbs. (57 kg)	Purifier Horizontal Clean Bench

*Enclosures supported by mobile Base Stands will require special modifications to incorporate flexible ducting, wiring and plumbing when applied.

Accessory Shelves

Glacier white, epoxy-coated, 16 gauge steel. Designed to hook on the lower rear horizontal cross member of the Telescoping Base Stand. Provides storage space for wipes, cleaning agents, gloves and other miscellaneous laboratory apparatus. One each.




Nominal Width	Catalog #	Shelf Width x Depth	Shipping Weight	For Use With:
2-Ft	3811100	20.4" x 9.9"	9 lbs. (4 kg)	3746700, 3746701
3-Ft	3811101	32.4" x 9.9"	12 lbs. (5 kg)	3746710, 3746711
4-Ft	3811102	44.4" x 9.9"	15 lbs. (7 kg)	3747002, 3746702, 3747012, 3746712
6-Ft	3811104	68.4" x 9.9"	23 lbs. (10 kg)	3746704, 3746714

ELECTRIC & MANUAL HYDRAULIC LIFT BASE STANDS

These stands have a built-in hydraulic lift so that height may be adjusted at any time. They allow you to position the stand and the safety enclosure or clean bench it supports to accommodate persons of different heights and wheelchair-bound users, to transport equipment through doorways, and for easier service access.

Standard Features

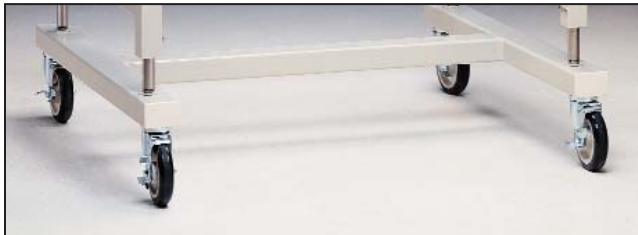- Support Purifier Class I Safety Enclosures, Purifier HEPA Filtered Enclosures, Purifier Horizontal Clean Benches, Purifier Vertical Clean Benches and Purifier Filtered PCR Enclosures.
- Support loads up to 1000 pounds.
- ADA-compliant.
- Durable 1.75" tubular, epoxy-coated steel frame with rear cross-bar supports. Sub-frame is 2.0" x 3.0".
- Height is infinitely adjustable from 25.5" to 33.5" high.
- Include four powder-coated, die-cast feet with levelers.
- Include front-accessible up/down switch (Electric Lift models)
- Include front-accessible hand crank. Turning the crank clockwise raises the stand; counterclockwise lowers it. The handle may be disengaged and folded under, out of the way, when not in use. (Manual Lift models)




Caster Kit 3784000

Casters add mobility to the Hydraulic Lift Base Stand. Kit includes four 5" diameter, non-marking polyurethane casters with toe locks. Load rating for each caster is 250 pounds. Casters have bearings for quiet operation. Casters replace the fixed feet that are included with the stand. **Installation is required.** Not for use with stands that support hard-ducted enclosures. (4.5 kg)




Manual Hydraulic Lift Base Stand 3780304.

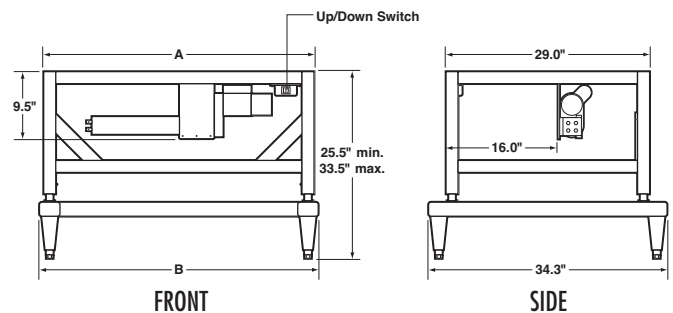Electric Lift



Manual Lift

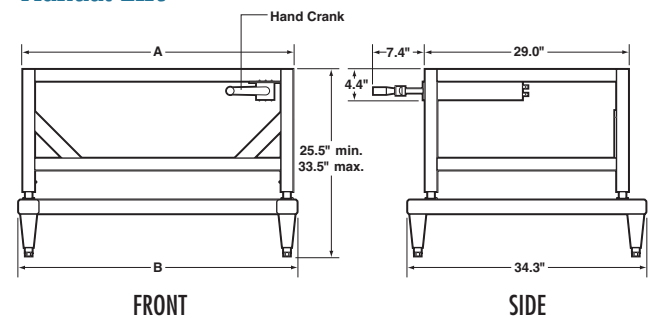

Catalog Numbers	Nominal Width	A	B
3780300, 3780303, 3780400	3-Ft	36.0"	37.2"
3780301, 3780304, 3780401	4-Ft	48.0"	49.2"
3780302, 3780305, 3780402	6-Ft	72.0"	73.2"

ELECTRIC & MANUAL HYDRAULIC LIFT BASE STANDS

Electric Hydraulic Lift Base Stands

Nominal Width	Catalog #	Overall Width x Depth	Switch Location	Electrical Requirements	Shipping Weight	For Use With:
3-Ft	3780300	36.0" w x 29.0" d	Right Side	115 volts, 60 Hz, 2.4 amps	150 lbs. (68 kg)	Purifier Class I Enclosure, Purifier HEPA Filtered Enclosure, Purifier Horizontal Clean Bench, Purifier Vertical Clean Bench, Purifier Filtered PCR Enclosure
3-Ft	3780303*	36.0" w x 29.0" d	Right Side	230 volts, 50 Hz, 0.8 amp	150 lbs. (68 kg)	Purifier Class I Enclosure, Purifier HEPA Filtered Enclosure, Purifier Horizontal Clean Bench, Purifier Vertical Clean Bench, Purifier Filtered PCR Enclosure
4-Ft	3780301	48.0" w x 29.0" d	Right Side	115 volts, 60 Hz, 2.4 amps	180 lbs. (82 kg)	Purifier Class I Enclosure, Purifier HEPA Filtered Enclosure, Purifier Horizontal Clean Bench, Purifier Vertical Clean Bench, Purifier Filtered PCR Enclosure
4-Ft	3780304*	48.0" w x 29.0" d	Right Side	230 volts, 50 Hz, 0.8 amp	180 lbs. (82 kg)	Purifier Class I Enclosure, Purifier HEPA Filtered Enclosure, Purifier Horizontal Clean Bench, Purifier Vertical Clean Bench, Purifier Filtered PCR Enclosure
6-Ft	3780302	72.0" w x 29.0" d	Right Side	115 volts, 60 Hz, 2.4 amps	220 lbs. (100 kg)	Purifier Horizontal Clean Bench
6-Ft	3780305*	72.0" w x 29.0" d	Right Side	230 volts, 50 Hz, 0.8 amp	220 lbs. (100 kg)	Purifier Horizontal Clean Bench

Manual Hydraulic Lift Base Stands

Nominal Width	Catalog #	Overall Width x Depth	Handle Location	Shipping Weight	For Use With:
3-Ft	3780400	36.0" w x 29.0" d	Right Side	150 lbs. (68 kg)	Purifier Class I Enclosure, Purifier HEPA Filtered Enclosure, Purifier Horizontal Clean Bench, Purifier Vertical Clean Bench, Purifier Filtered PCR Enclosure
4-Ft	3780401	48.0" w x 29.0" d	Right Side	180 lbs. (82 kg)	Purifier Class I Enclosure, Purifier HEPA Filtered Enclosure, Purifier Horizontal Clean Bench, Purifier Vertical Clean Bench, Purifier Filtered PCR Enclosure
6-Ft	3780402	72.0" w x 29.0" d	Right Side	220 lbs. (100 kg)	Purifier Horizontal Clean Bench

* International electrical configuration

ACCESSORIES

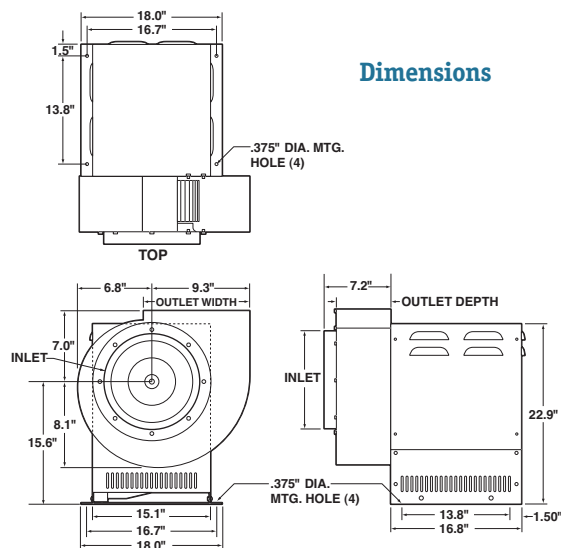


Coated Steel Blowers for Canopy-Ducted Purifier Class I and HEPA Filtered Enclosures

A remote blower or house exhaust is required when canopy-ducting the Purifier Class I Safety Enclosure or Purifier HEPA Filtered Enclosure.

Phenolic-coated impeller and housing, dry powder epoxy-coated steel base and weather cover, belt-drive

motor with adjustable sheave and self-adjusting gravity belt tightener, flow-through vent holes and louvers, lubricated pillow block bearings, and variable exhaust orientation (vertical discharge is recommended). Blower inlet is 10.87" OD and is sized to accept 10" nominal diameter PVC duct. Blower outlet is 5.5" d x 10.0" w. Contact Labconco for ordering information on Remote Blowers with other equivalent duct lengths or diameters or other electrical configurations.



Dimensions

Catalog #	For use with:	Electrical Requirements	HP	.12* CFM@RPM	.25* CFM@RPM	.38* CFM@RPM	.50* CFM@RPM	.62* CFM@RPM	.75* CFM@RPM	Shipping Weight
7068000	Canopy-ducted 2-Ft & 3-Ft Purifier Class I & HEPA Filtered Enclosures with up to 100 equivalent feet of 6" dia. duct	115 volts, 60 Hz, 4.0 amps	1/6	370 @ 530 500 @ 650	300 @ 635 540 @ 800	308 @ 760 410 @ 800	—	—	—	92 lbs. (41.7 kg)
7068200	Canopy-Ducted 4-Ft Purifier Class I & HEPA Filtered Enclosures with up to 100 equivalent feet of 6" dia. duct	115 volts, 60 Hz, 4.4 amps	1/4	—	540 @ 800 720 @ 950	410 @ 800 760 @ 1050	350 @ 870 710 @ 1060	390 @ 970 620 @ 1060	430 @ 1060	92 lbs. (41.7 kg)

* Static pressure in inches H₂O

ERGONOMIC CHAIR




The Ergonomic Chair provides comfortable seating and may be used throughout the laboratory.

Standard Features


- 6-way articulating seat and back control for personalized adjustment.
- Pneumatic mechanism seat height adjustment from 18.25" to 25.75".
- Five-leg black reinforced composite base with 2" diameter ball-bearing casters.
- Aluminum support ring.
- Removable armrests.
- Black vinyl upholstery.

Catalog #	Description	Shipping Weight
3744000	Ergonomic Chair with Armrests	35 lbs. (15.9 kg)

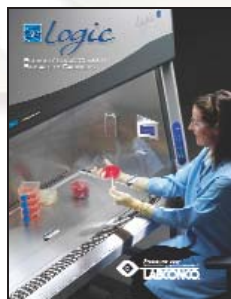*Contact Labconco for more information about these
quality products for your laboratory.*




Biological
Safety Cabinets

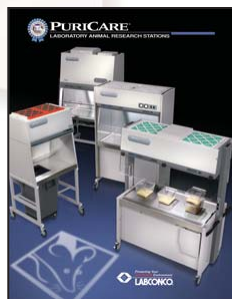


Laboratory Animal
Research Stations



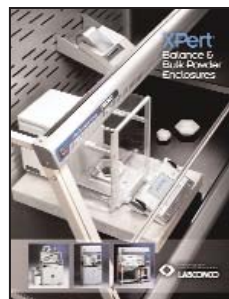Fume Hoods &
Carbon-Filtered
Enclosures



Balance & Bulk Powder
Enclosures

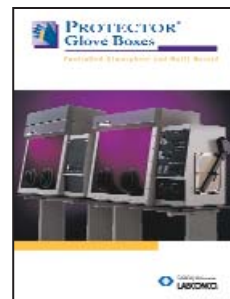

Glove Boxes



Glassware
Washers



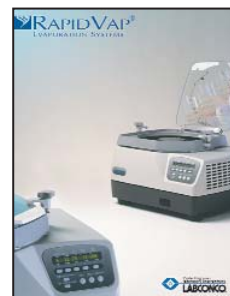
Water Purification
Systems



Freeze Dry
Systems



Centrifugal Concentrators
& Cold Traps




Multiple Sample
Evaporation Systems



Agricultural Chemistry
Products



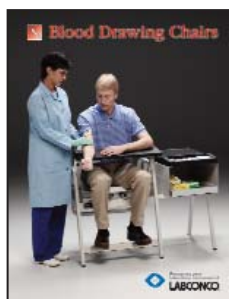Laboratory
Carts & Benches



Blood Drawing
Chairs



Digital
Chloridometers



Vacuum
Desiccator



Labconco Corporation

8811 Prospect Avenue
Kansas City, MO 64132-2696
816-333-8811 or 800-821-5525
FAX: 816-363-0130
E-MAIL: labconco@labconco.com
HOME PAGE: www.labconco.com



© 2007 by Labconco Corporation
Printed in the U.S.A.
Product design subject to change without notice

2-12-9/07-GC-10M-R1